


KDHE Microbiology Lab

Packaging and Shipping Guide for Water Samples Collected for Microbiological Analysis

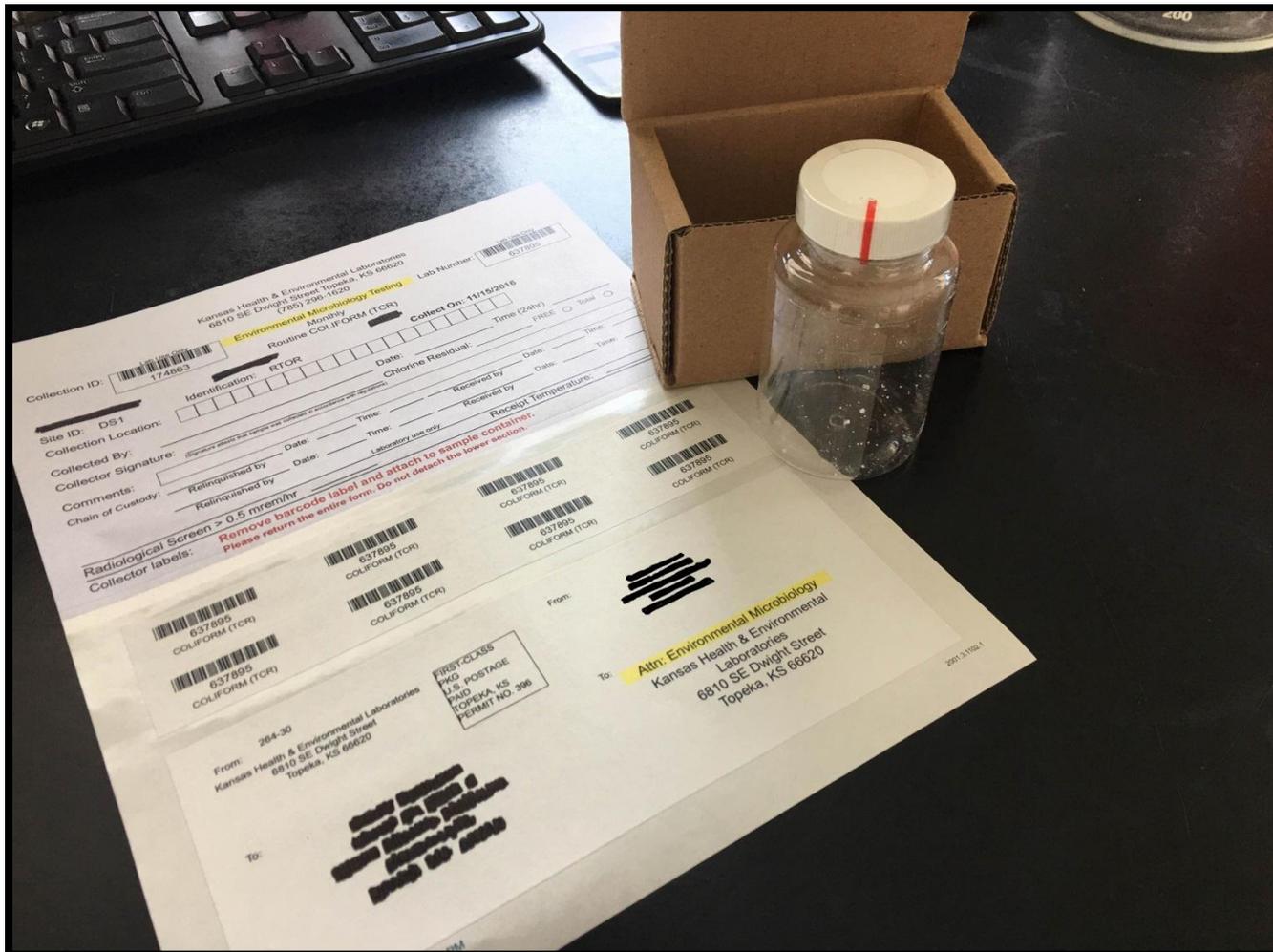


Collection instructions are included with your sample bottle kit and can be found at www.kdheks.gov/envmicro

The shipping box **should be** reused to send sample back for analysis



Collection kits contain the following items:

1. Sterile Idexx collection bottle
2. Collection form with collection site, information, and barcodes
3. Return address label
4. Yellow collection information sheet
5. Instructions for specific sample type (e.g. Replacement, Repeat, Triggered Ground Water)

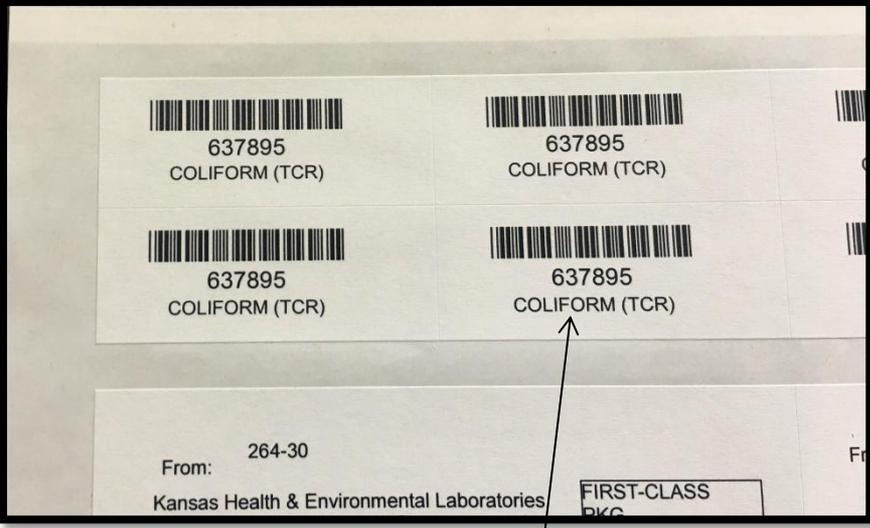

3. Make sure to remove a barcode label..

And attach it to the sample bottle vertically

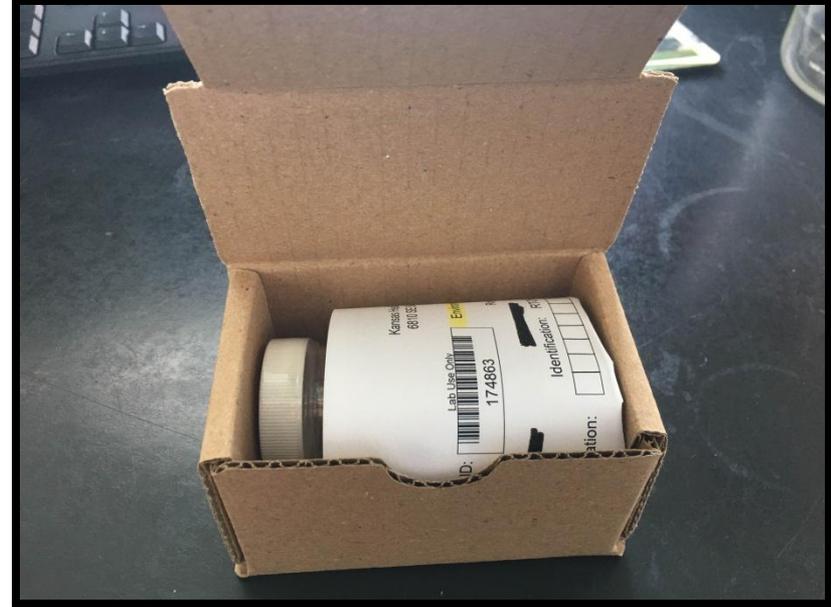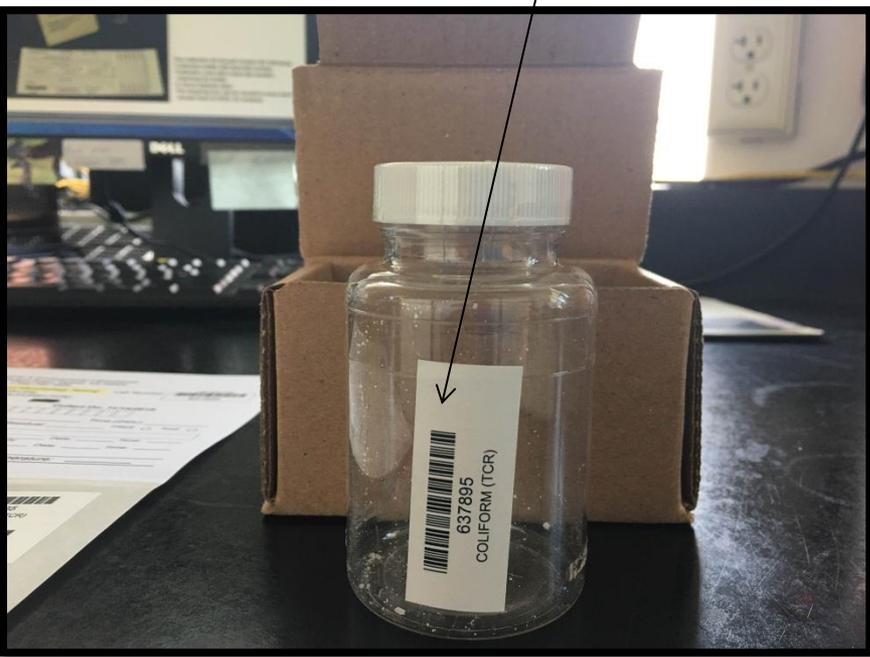

4. Wrap the completed form around the bottle and place in the box

5. Attach return label to box and mail to KHEL. All done!

